

Quick Guide to FluorPen Operation

1. To Switch on: Hold the Set key for one second.
2. To Scroll Through a Menu: Use the Menu key.
3. To Select a Menu Option: Use the Set key.
4. To Get Into a Submenu: Scroll through the main menu and find your option based on the cursor (>). Use the Set key to make your selection and to get into a submenu.
5. To Scroll Through a Submenu: Use the Menu key.
6. To Return From a Submenu: Scroll through the submenu up to the >Return option. Use the Set key to make your selection and to return to the main menu.
7. To Switch off: Scroll to the >Turn Off command in the main menu and click the Set key. (The FluorPen switches off automatically after five minutes without a command).

Detailed instructions with step-by-step guiding through all the FluorPen functions are in your owner's Manual.

Quick Guide to FluorPen Operation

Main Menu:

Measure
Data
Setting
Turn Off

Main Menu with SubMenus:

Measure
- *Ft*
- *QY*
- *Return*
Data
- *Browse*
- *Erase*
- *Return*
Setting
- *f-pulse*
- *F-pulse*
- *Sound*
- *Return*
Turn Off